
Cell Freezing Container

BS-05-CFC For 12 x 5mL Cryogenic Vials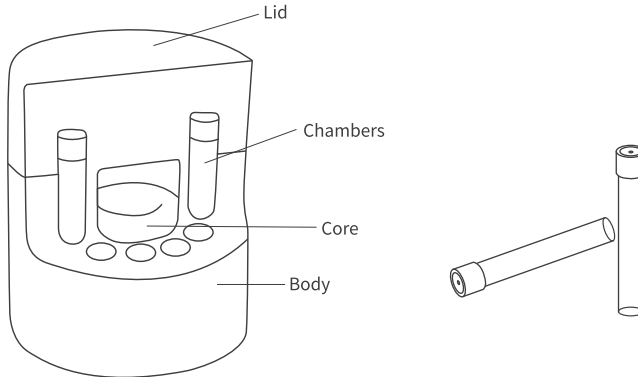

Biosharp Cell Freezing Container, can provide alcohol-free freezing at the rate of $-1^{\circ}\text{C}/\text{minute}$ that is ideal for cryopreservation of most cells and cell lines with a -80°C freezer. Cell Freezing Container is made of cross-linked polyethylene foam with a solid state core (metal ring). The foam is non-absorbent and won't affect nearby frozen samples. Freezing profiles are consistent and reproducible.

Features:

- Ease to use
- Alcohol and fluid-free freezing
- Lower cost of use than alcohol-based devices
- High cell recovery and cell viability
- Reproducibility

Instruction:

1. Place sample vials containing cell suspension in each well of the Cell Freezing Container under room temperature. Fully seat the lid on Cell Freezing Container.
2. Place Cell Freezing Container upright into a -80°C freezer. Freeze for minimum four hours to achieve standard cryopreservation effect.
3. Take out Cell Freezing Container and transfer vials immediately to Liquid Nitrogen or for other following applications.
4. Place Cell Freezing Container in room temperature for next freezing until the foam and core are fully return to room temperature.

Cautions:

1. Ensure that there is at least 2.5cm of free space clearance around Cell Freezing Container when placing in the -80°C freezer.
2. Make sure the core is seated in the bottom of the central cavity, and lid is fully seated on the container before placing in the freezer.
3. Maximum temperature exposure: 60°C . Avoid prolonged exposure to UV light sources.
4. It is strongly recommended that all frozen cell cultures be checked for viability before the stock culture is terminated.

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